

Morphologi G3 – Malvern Panalytical

INTRODUCTION

The Lumetics LINK™ software platform scans network locations for new measurement data files, copies data directly to a centralized database, and provides a powerful user interface for rapid multi-measurement multi-technique data aggregation, visualization, analysis, and reporting. LINK employs a client/server-based architecture where the LINK server hardware is provided by the end user and resides on the end user's network. The LINK client is a portable web-based application that may be placed on any computer with network connectivity to the LINK server. For successful import, the LINK webserver requires read access to the folders where user data resides.

The Morphologi G3 provides an advanced particle characterization tool for the measurement of particle size and particle shape from 0.5 microns to several millimeters.

DETAILS

LINK requires the exported.TXT and particles.TXT data files. The BMP particle image files are optional. Data files must be exported from the Morphologi G3 software using the Lumetics G3 Export Template.EDF, which is available in the customer portal on the Lumetics Website (http://lumetics.com/customer_portal/).

The following raw curve data may be imported, in addition to all available instrument/analysis settings and parameters calculated by the instrument software:

- Particle count or concentration vs. morphology

Particle count/concentration for each supported morphological parameter occurs at a pre-defined bin spacing and measurement range as per the table in the Users manual (section 4.29).

LINK requires the data to be in the following format:

- One folder per measurement
- Metadata is exported in the *exported.txt* file, with the Record # as the first column
- Particle information is found in the *particles.txt* file
- Thumbnails are exported in the same folder as the *exported.txt* and *particles.txt* folder.

The following is the default Morphologi G3 specific export method:

1. Create Folder for Exported Data
 - In Windows, create a folder for each of the exported Morphologi G3 files.
 - NOTE: Each exported file must be in its' own folder.
2. Copy LINK Export Template (NOTE: Perform this step once)
 - Contact Lumetics for the *LINK G3 Export Template (.edf)* file.
 - Copy this file to the Documents folder containing the Morphologi G3 Export Templates.
 - By default, this will be: C:\Users\Public\Documents\Malvern Instruments\Morphologi\Export Templates
 - NOTE: the location may vary depending on installation preferences (e.g. it may be in a specific user's folder).
3. Export Results
 - In the Morphologi G3 Records tab, select a single measurement for export.
 - Select File -> Export -> Export Result... and configure menu as shown below.
 - Browse for the location of the data folder (see Step 1 above). The filename MUST be 'exported.txt'
 - In the Parameters tab, select Use export template parameters and select the Lumetics G3 Export Template. If you do not see this template, return to Step 2.

4. Export Particles

- From the Morphologi G3 main screen, first ensure the record is still selected and then select File -> Export -> Export Particles...
- Browse to folder from Step 1, name the file particles.txt (MANDATORY) and select other options as shown below in File tab. Leave all other tabs as Default.


5. Export Images

- From the main screen of the Morphologi G3 software select File -> Export -> Export Images.... Select the data folder from Step 1. Click OK.
- Note: it may take a few minutes to export the images.

EXAMPLES


Included below are sample dashboards from Morphologi G3 measurement files:

1. Radar Chart plotting multiple particle characterization parameters with limit lines/acceptance mask

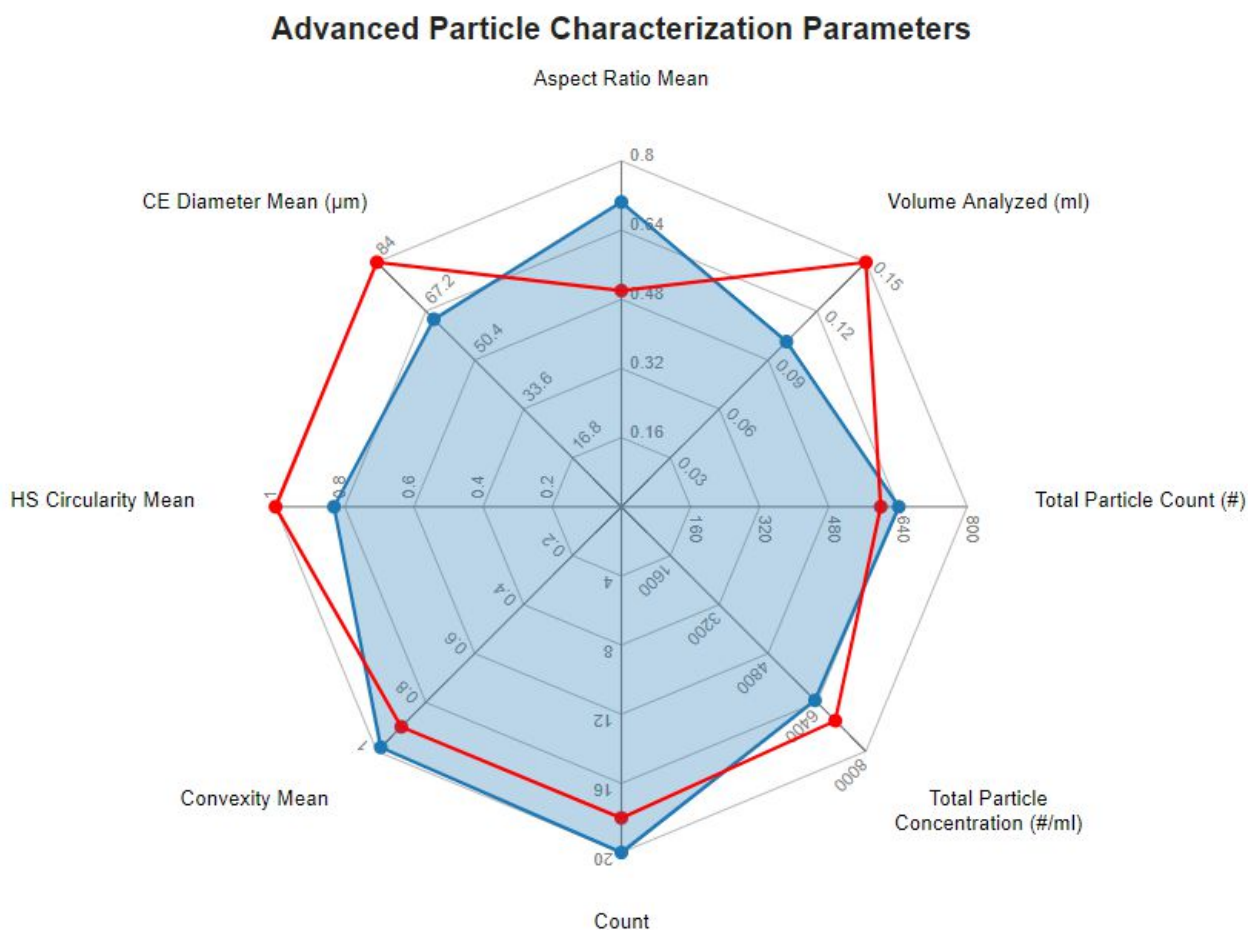

Morphologi G3, Total Population

2. Histogram plotting binned raw data for Particle Concentration vs. Diameter

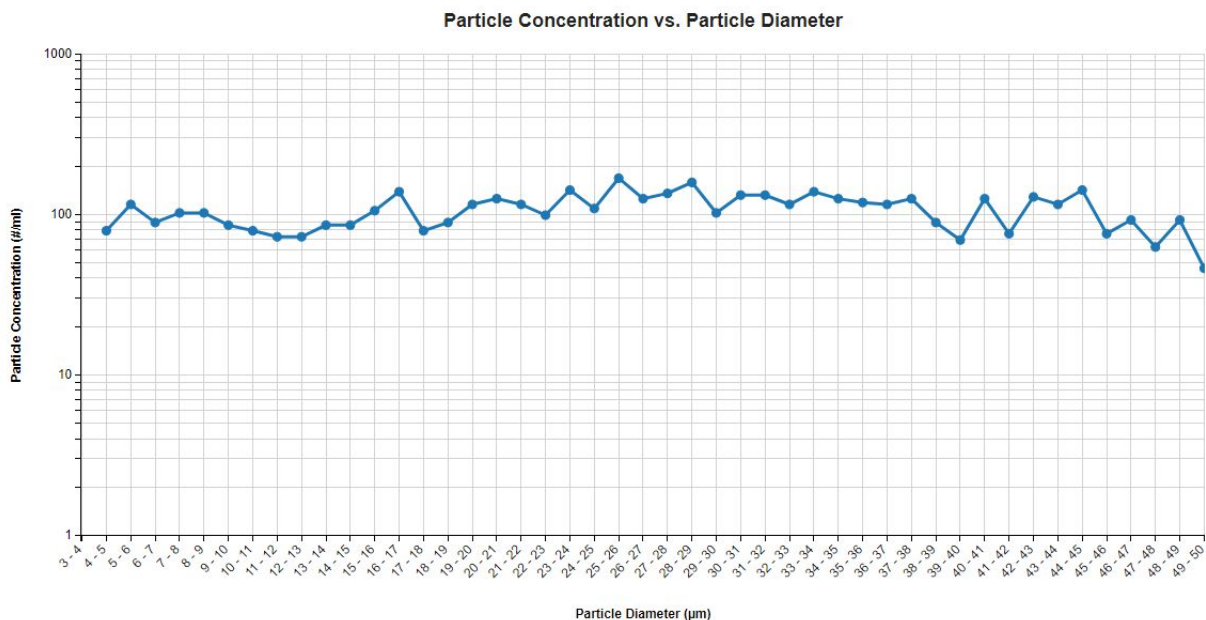

3. Tabular Summary examples

Measurement Summary Table – Measurement Results

Sample Name	Volume Analyzed (ml)	Volume (%) (Particle Diam. (µm) >=1) - AVG	Particle Count (#) (Particle Diam. (µm) >=1) - AVG	Total Particle Concentration (#/ml) - AVG	Total Particle Count (#) - AVG	≥1µm AVG	2-10µm - AVG	≥10µm AVG	HS Circularity Mean - AVG	Count - AVG	Convexity Mean - AVG	CE Diameter Mean (µm) - AVG	Aspect Ratio Mean - AVG
Sample A	0	100	1	9.87	1	10	0	0	0.84	15	0.93	62.00	0.80
Sample B	0	100	1728	17058.24	1728	17,058	1,718	15,341	0.82	22	0.89	60.00	0.50
Sample C	0	100	196	1934.85	196	1,935	0	1,935	0.83	20	0.98	64.43	0.71

Measurement Summary Table – Instrument Settings

Sample Name	Analysis Date	Version	Source ID - AVG	S/W Version	Record # - AVG	Instrument S/N	Image DB	Date
Sample A	2017-01-31 15:4...	8	10	8	10	MAL1147555	Training_10(1).img	31 January 2017 1...
Sample B	2011-05-25 11:2...	7.40.029	2	7.40.029	3	MAL1009013	Example Data_2(1...	May 25, 2011 11:...
Sample C	2012-08-29 08:0...	7.40.029	5		8	MAL1009015		August 29, 2012 8...

DASHBOARD DOWNLOAD

Included below is a link to downloadable dashboards for Morphologi G3 measurement files:

<http://lumetics.com/dashboards/ MorphologiG3/MorphologiG3.zip>

CONTACT LUMETICS

For direct assistance, please contact Lumetics LINK™ Support:

E-mail: support@lumetics.com

Phone: 1.613.614.874

Website: <http://lumetics.com/>

